Increasing Efficiency and Pesticide Recovery from the QuEChERS Approach for Fruit and Vegetable Products Using Ceramic Homogenizers

Joan Stevens, Laure Bertin, and Melanie Rothermich, Agilent Technologies, 2850 Centerville Road, Wilmington, DE, 19808, USA

Introduction

QuEChERS

- > Quick, Easy, Cheap, Effective, Robust and Safe
- Developed by the US FDA and EU Food Regulatory Agencies
- Procedure was validated in 2003, "toddler stage"
- > Extraction and analysis of pesticides in food product
- > Advancements in the analysis of other components
- > PCB, PAH, Antibiotics, Pharmaceuticals, Toxins
- \succ Other food-stuffs; meat fish, grains, nuts, juices, oils
- \succ Other matrices; soil, wine, biologicals

➢ QuEChERS

- Basic procedure
- Concerns and variance with procedure

Homogenizers in QuEChERS extraction

- Description and characterization
- > Implementation
- Data and results

Experimental

QuEChERS Procedure

> 3 step process: extraction, dispersive-SPE, analysis



 \geq 2nd Dispersive-SPE





Concerns and Variance with Procedure

- > 1st Step Extraction
- > After addition of organic to sample, add extraction salts
- > Requires vigorous shaking, 1 min or greater
- > Variations is time and force shaking affect recovery
- \succ Variation from beginning to end of extraction, 100+ samples
- > Variations between technicians between labs
- > Affect recovery of pesticides

Homogenizers in QuEChERS Procedure

- Homogenizers are inert ceramic pieces
- QuEChERS
- Break up salt agglomerates

Ceramic Homogenizers (CH)

- \blacktriangleright Available in 3 sizes; 50 mL, 15 mL, and 2 mL
- ➢ Place 2 CH per tube
- Process samples the same
- shaking capabilities

- ► AOAC method 2007.01 \succ 6 g MgSO₄, 1.5 g NaAcetate
- \geq EN method 15662



EPRW 2010

Experimental

Results and Discussion

> Aggressively force the partitioning between the water and organic layers in

Homogenizers

Increase grinding/pulverizing of homogenized matrix



> Weight and angled sides of the ceramic homogenizers increase



QuEChERS Extraction: With and Without Ceramic

> Determination of inertness of CH in the presence of matrix and pesticides

> Study to extract samples with and without CH by manual shaking

 \rightarrow 4 g MgSO₄, 1 g NaCl, 1 g NaCitrate, 0.5 g disodium citrate sesquihydrate

Results and Discussion



Graph 1: Pesticides used in QuEChERS study with and without ceramic homogenizers

Procedure after Addition of Ceramic Homogenizers

> 1st Step Extraction

With Ceramic Homogenizers

- > After addition of organic to sample, add extraction salts
- ✓ Requires vigorous shaking, 1 min or greater
- ✓ Variations is time and force shaking affect recovery
- ✓ Variation from beginning to end of extraction, 100+ samples
- ✓ Variations between technicians between labs ✓ Affect recovery of pesticides
- "SHAKING"

- > Maintain acceptable recovery and RSD for all pesticides, classes \succ Setup a timed extraction study
- \succ Add the extraction salts

Average Recovery of Several Classes of Pesticides within 20 Second Shaking **Extraction and Ceramic Homogenizers**



Graph 2: Pesticides used in study: Acephate, Carbaryl, Carbendazim, Cyprodinil, Imidacloprid, Imidazalil, Methamidophos, Penconazole, Propoxur, Pymetrozine, Thiabendazole, Thiophanate-methyl, Ethoprophos, Kresoxim-methyl ; Apple matrix

Conclusions

- \succ QuEChERS procedure is very efficient "just enough" sample preparation
- Ceramic homogenizers addresses many of the "shaking" issues
- > Inert, improves recoveries and RSDs for the pesticides and other compounds
- > Disposable, can be used manually or with mechanical shakers
- Drastically decreases "shaking" time by 70%
- Improve sample throughput

Agilent Technologies

- AOAC Method with Ceramic EN Method with Ceramic Homogenize
- AOAC Method without Ceramic EN Method without ceramic

QuEChERS Extraction: Require Shaking for 1 min

Investigate the ceramic homogenizers

- \succ Could they reduce the time required for shaking
- > Shake for 10 seconds, adding 10 sec for each additional set
- > 10, 20, 30, 40, 50, 60 sec